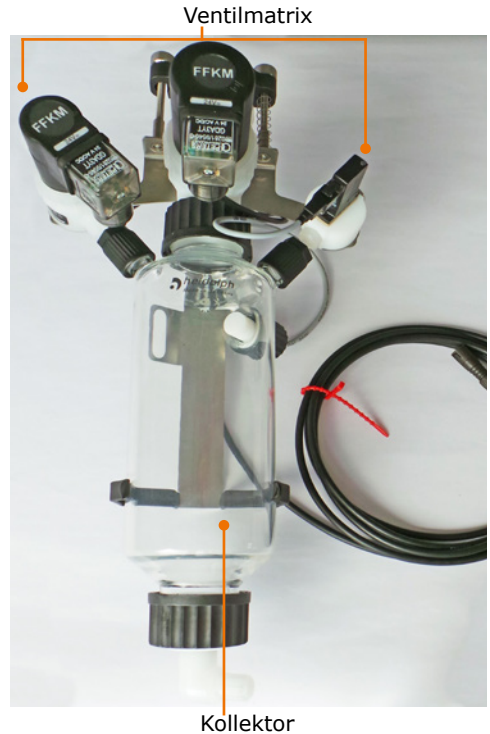
➤ Ventilmatrix und Kollektor

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Best. Nr. 569-01000-00

- 1x Kollektor
- 1x Ventilmatrix

Ventilmatrix

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
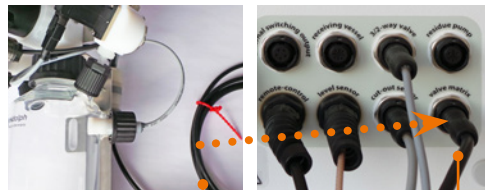
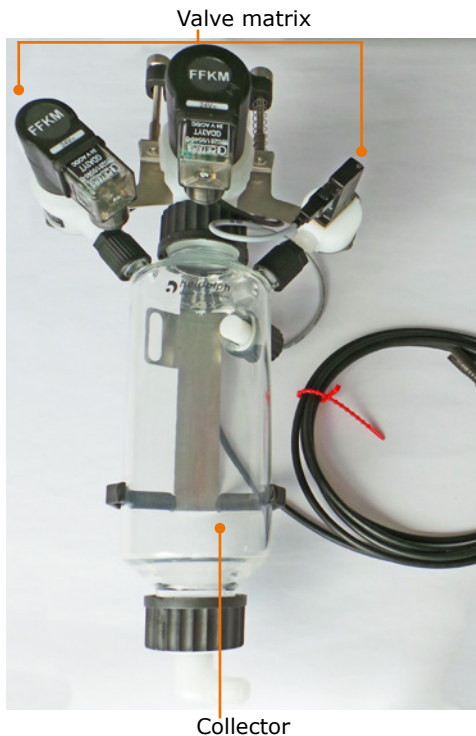
- 1x Collector
- 1x Valve matrix

Valve matrix



The valve matrix on the rotary evaporator controls evacuation and ventilation of the system as well as emptying the condensate collector.

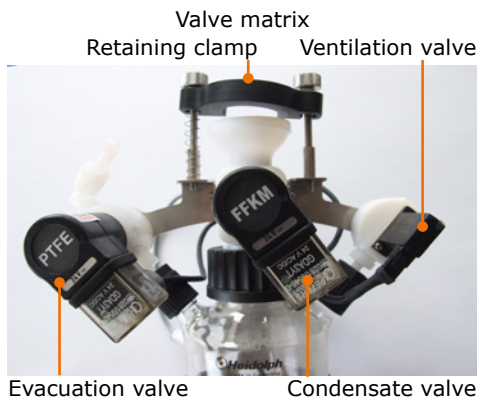- If necessary remove the receiving flask from your rotary evaporator.
- Open the retaining clamp at the valve matrix.

Connecting valve matrix to Hei-VAP Industrial

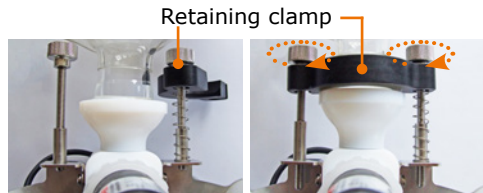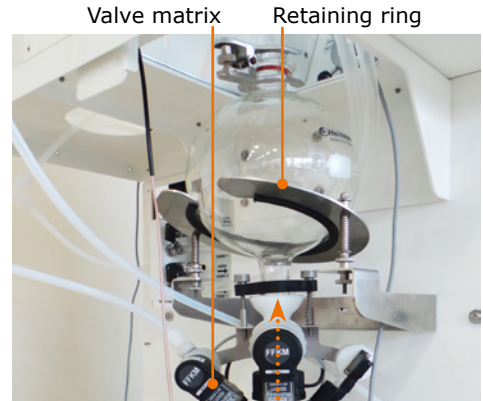- If necessary remove the cassette system including receiving flask from the Hei-VAP Industrial.
- Mount the holdig system for the buffer flask instead.
- Press the retaining ring downwards against the spring force and place the buffer flask on top of the retaining ring.
- Release the spring force carefully until the upper flange touches the glass condenser.
- Connect buffer flask and glass condenser using the ground joint clamp.
- Guide the condensate collector from below to the buffer flask.
- Close the retaining clamp using the two knurled screws on the right and left.
- Tighten the two knurled screws evenly.

Connecting valve matrix to control box

All cables of the valve matrix are firmly pre-assembled. They are joined at the lower outlet to form a control cable.

- Connect the valve matrix control cable to the "valve matrix" connector on the back of the control box.

Valve matrix control cable

Valve matrix connector