Safe  
Heating and Mixing

## Safety Guide

Magnetrührer  
Magnetic stirrers

## Index

German.....	3	
English.....	10	
Bulgarian .....	17	
Czech .....	24	
Danish .....	31	
Greek .....	38	
Español .....	45	
Eesti .....	52	
Suomi .....	59	
French .....	66	
Hrvatski .....	73	
Magyar .....	80	
Italian.....	87	
Lithuanian .....	94	
Latviešu .....	101	
Nederlands .....	108	
Polski.....	115	
Português.....	122	

Romanian.....	129	
Slovensky.....	136	
Slovenian.....	143	
Svenska.....	150	

## Zu diesem Dokument

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## About this document

This document contains a summary of all safety instructions that must be observed when handling magnetic stirrers from Heidolph instruments GmbH & Co. KG.

This document in no way replaces the operating instructions for the product!

For more detailed information, please refer to the original instructions (German language) or the translation of the original instructions available on demand.

## Typographic conventions

Standardized symbols, highlighting elements, and signal words are used in this document to identify warnings, cautions, important information, and special text contents.

Symbol	Signal word / explanatory note
	Warning symbols in combination with a signal word indicate dangers: <b>DANGER</b> Indicates a hazardous situation which, if not avoided, will result in death or serious injury.
	<b>WARNING</b> Indicates a possible hazardous situation which, if not avoided, could result in death or serious injury.
	<b>CAUTION</b> Indicates a hazardous situation which, if not avoided, could result in minor or moderate injury or material/environmental damage.
	Mandatory signs are used to indicate important information regarding the product handling. This information is used to ensure operational safety and to maintain the value of the product.

## Copyright protection

This publication is protected by copyright and intended for internal use by the purchaser of the product only.

No part of this publication may be transmitted or reproduced in any form, by any means, without the prior written consent of the copyright owner Heidolph Instruments GmbH & Co. KG. Any violation is subject to compensation for damages.

## Intended use

Magnetic stirrers from Heidolph instruments GmbH & Co. KG are specifically designed for the following laboratory tasks:

- Heating
- Mixing
- Stirring
- Titration

Any other use of the device described in this manual is not considered as intended!

Due to its design, the device in its delivery condition may only be used in analytical processes or in laboratory-like conditions in the food, cosmetics, and pharmaceutical industries as well as other comparable industries that manufacture products intended for consumption by humans or animals, or for use on humans or animals.

## General safety information

- Before commissioning and using the device, familiarize yourself with all the safety regulations and guidelines for occupational safety applicable at the place of use and observe them at all times.
- Only operate the device if it is in perfect technical condition. In particular, ensure that there is no visible damage on the device itself and, where necessary, on connected devices or the supply connections.
- If there is missing or misleading information on the device or on occupational safety, contact the responsible safety specialist or our technical service.
- Only use the device in accordance with the regulations for intended use.

## Electrical safety

- Ensure that the voltage indicated on the rating plate matches the supply voltage of the country in which the device is being used.
- Ensure that the mains socket-outlet is protected by means of a residual-current device (RCD).
- Always use the supplied power supply 3-pin cord provided with the device (phase, neutral, ground).
- Prior to use, check that the device and the power supply cord are free of visible damage.
- Have repairs and/or maintenance work on the device carried out exclusively by an authorized and skilled electrician or by the technical service department of Heidolph Instruments.
- Always switch the device OFF and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

## Operational safety

- Operate the device under a closed ventilated fume hood when working with potentially hazardous substances (see EN 14175 and DIN 12924).
- Do not make any unauthorized changes or modifications to the device!
- Only use genuine spare parts and accessories, or those expressly approved by the manufacturer!
- Rectify malfunctions or faults on the device immediately.
- Shut down the device and disconnect it from the power supply if it is not possible to eliminate the malfunction or rectify the fault immediately.

- Observe all other applicable regulations such as laboratory and workplace guidelines, recognized safety technology rules and special local regulations.

## Personal protective equipment (PPE)

The operator must determine and provide the necessary PPE, depending on the respective application and the media and chemicals used.

The corresponding instruction of the personnel is solely within the operator's responsibility.

## Environmental protection

When processing environmentally hazardous substances, take appropriate measures to avoid risks to the environment.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the responsible personnel is the sole responsibility of the operator!

## Biohazard

When processing biohazardous substances, take appropriate measures to prevent hazards to persons and the environment, including:

- Instruction of the personnel regarding the necessary safety measures.
- Provision of personal protective equipment (PPE) and instruction of the personnel in its use.
- Marking of the device with a biohazard warning symbol.

The evaluation of corresponding measures such as the marking of a hazardous area, their implementation, and the training of the responsible personnel is the sole responsibility of the operator!

## Other regulations

In addition to the notes and instructions in this document, observe all other applicable regulations such as laboratory and workplace guidelines, hazardous substances ordinances, recognized rules of safety engineering and occupational medicine as well as particular local regulations!

## Work safety

### General instructions

- Always use the prescribed personal protective equipment (PPE) such as protective clothing, safety goggles, protective gloves, safety shoes, etc.
- Persons with pacemakers must maintain a minimum distance of 10 cm from the device to prevent possible interference.
- Do not operate any other devices in the immediate vicinity of the device ...
  - which can generate electromagnetic fields in the frequency range between  $9 \times 10^3$  Hz to  $3 \times 10^{11}$  Hz,
  - which generate emission or radiation sources in the frequency range  $3 \times 10^{11}$  Hz to  $3 \times 10^{15}$  Hz (in the optical spectral range wavelengths from 1,000  $\mu\text{m}$  to 0,1  $\mu\text{m}$ ),
  - which generate ultrasonic or ionizing waves.

- Do not process substances that could release energy in an uncontrolled manner (e.g. self-ignition).
- Do not process substances in which the energy input through mixing poses a danger.
- Do not work with an open flame in the immediate vicinity of the device (risk of explosion).
- Wipe off any fluid that may have spilled on the device immediately.
- Always switch off the device when it is not in use.

## Assembly

### CAUTION

#### Risk Of Slipping!

The vibration during operation can cause the device to move over the mounting surface and/or fall to the ground.

- Observe the instructions for the correct positioning of the device!

## Operation

### CAUTION

#### Risk of injury, risk of damage to property due to tilting/falling of the device

High shaking frequencies in combination with a large load and/or a tall attachment assembly on the shaking platform may cause the device to swing up and fall!

- Only gradually increase the shaking frequency to the required level under the above conditions, paying attention to the stability of the structure.
- Lower the shaking frequency or reduce the total load on the shaking platform when the device starts to move during operation.

An uneven distribution of the load on the shaking platform may cause the device to swing up and fall!

- Always ensure that the vessels are evenly distributed on the shaking platform.
- Pay particular attention to an even distribution of the weight load when vessels of different sizes and/or differently filled vessels are placed on the shaking platform at the same time!

#### Risk of injury, risk of damage to property from falling vessels

At high shaking frequencies, there is a risk that improperly fastened vessels will fall off the shaking platform.

- Before switching on the device, make sure that all vessels are properly fixed on the shaking platform.
- Contaminated surfaces should be cleaned immediately.

## Automatic restart

---

### WARNING

#### Risk of burns, risk of injury, risk of damage to property from falling vessels and leaking chemicals

When the [automatic restart] function is activated, the heating plate automatically heats up to the last set operating temperature when the power is restored and the rotation speed will be augmented to the last set value.

- Ensure that all vessels are properly positioned on the heating plate before the power is restored.
  - In the event of a power failure, do not place any objects on the heating plate.
- 

## Heating plate, heating function

---

### WARNING

#### Risk of burn

The heating plate is designed for a maximum operating temperature of 300°C. Contact with surfaces above 50 °C can lead to serious injury.

- Always pay attention to the heating temperature indicator of the device.
  - Avoid direct skin contact with the heating plate during operation and also observe the residual heat indicator after switching off!
  - Do not place heat-sensitive objects on the heating plate.
  - Always use the necessary personal protective equipment (heat-resistant gloves, eye protection, safety clothing) to process samples from a temperature of 50 °C.
- 

## Residual heat indicator

---

### WARNING

#### Risk of burn

As long as the device is not switched off, the display shows the current temperature of the heating plate or the measured value from the external temperature sensor.

- Note that the sensor reading may differ significantly from the current surface temperature of the heating plate!
  - Avoid direct skin contact with the heating plate!
  - Do not place heat-sensitive objects on the heating plate.
-

## Mixing function, rotation speed

---

### **WARNING**

#### **Risk of injury, risk of damage to property caused by splashing fluids**

When using open vessels, there is a risk that fluid will spray out.

- Whenever possible, use closed vessels for processing corrosive, toxic or biohazardous substances and seal them safely.
  - Especially using open vessels, increase the rotation speed only gradually and observe the fluid movements.
  - Always use the necessary personal protective equipment (heat-resistant gloves, eye protection, safety clothing) to process samples from a temperature of 50 °C.
- 

## Operation with external temperature sensor/temperature control

---

### **WARNING**

#### **Risk of burn**

As long as the temperature sensor is not immersed in the sample, the temperature of the ambient air is displayed. The heating plate can therefore heat up to 300 °C unnoticed!

- Always immerse the temperature sensor in the sample before switching on the device or before calibrating!
  - Do not place heat-sensitive objects on the heating plate.
  - Always use the necessary personal protective equipment (heat-resistant gloves, eye protection, safety clothing) to process samples from a temperature of 50 °C.
-

## Device service

---

### **WARNING: Danger of electric shock**

Live components are installed inside the device.

When opening the device, there is a risk of touching live components.

→ Switch the device's main switch off and disconnect it from the power supply before carrying out maintenance work, cleaning, or repairs.

Penetrating liquid poses the danger of an electric shock.

→ When cleaning, avoid the penetration of liquids.

---

## General cleaning instructions

---

### **CAUTION: Damage to the device**

Improper cleaning can damage the surfaces of the device.

Penetrating liquid can damage the electronic components inside the device.

→ Clean the device's surfaces with a soft, lint-free and only slightly moistened cloth.

→ Never use any aggressive or abrasive cleaning agents or aids.

---

## Repairs

Repairs to the device may only be carried out by authorized skilled experts!

Unauthorized repairs during the warranty period will result in the loss of the warranty claim.

The owner is solely liable for damage caused by unauthorized repairs.

In case of repair contact an authorized dealer or our technical service.

## Maintenance

The device contains no user-serviceable components. If necessary, in the event of abnormal operating behavior such as excessive noise or heat generation, for example, contact our technical service.

## Dismantling

Observe the instructions contained in the assembly instructions supplied when dismantling.

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© Heidolph Instruments GmbH & Co. KG